



MicroVue Bone Rat PTH EIA Sample Diluent

QUIDEL

For **Research Use Only**. Not for use in diagnostic procedures.

A symbols glossary can be found at quidel.com/glossary

This reagent is intended for research use only in the determination of rat parathyroid hormone. It should be used to dilute serum or plasma samples with intact and/or fragment PTH concentrations exceeding the measurable range of the immunoassay. To obtain final results multiply the observed concentration by the sample dilution factor. [i.e. 180 pg/mL (observed) x 10 (dilution factor) = 1800 pg/mL].

This sample diluent should only be used when measuring samples in either the Immotopics Rat Intact PTH ELISA Kit #60-2500 or Rat Bioactive Intact PTH ELISA Kit #60-2700. It is not an appropriate sample diluent for use in any other kit.

PREPARATION AND STORAGE

Store this reagent at 2°C to 8°C upon receipt. One bottle containing 10 mL of a lyophilized, treated, stabilized, protein matrix with 0.1% Ciprofloxacin as preservative. Before use, allow to come to room temperature and reconstitute with 10 mL of deionized water. Allow the bottle to sit for approximately 20 minutes with occasional gentle swirling and inversion. Assure complete reconstitution before use.

After reconstitution the sample diluent may be stored at 2°C to 8°C for up to 3 days. If longer storage is necessary the diluent should be aliquoted in volumes to minimize freeze/thaw cycles and stored frozen at –20°C or below. The reagent is stable until the expiration date on the label.

WARNINGS AND PRECAUTIONS

- In case of contact with the sample diluent, wash thoroughly with water.
- Use Good Laboratory Practices.
- Wear suitable protective clothing, gloves, and eye/face protection when handling the contents of this kit.
- Dispose of containers and unused contents in accordance with Federal, State and Local regulatory requirements.
- Wash hands thoroughly after handling.
- For additional information on hazard symbols, safety, handling and disposal of the components within this kit, please refer to the Safety Data Sheet (SDS) located at quidel.com.

ASSISTANCE

To place an order or for technical support, please contact a Quidel Representative at 800.874.1517 (in the U.S.) or 858.552.1100 (outside the U.S.), Monday through Friday, from 8:00 a.m. to 5:00 p.m., Eastern Time. Orders may also be placed by fax at (740) 592-9820. For e-mail support contact customerservice@quidel.com or technicalsupport@quidel.com.

For services outside the U.S.A., please contact your local distributor. Additional information about Quidel, our products, and our distributors can be found on our website quidel.com.

REF

30-2531 MicroVue Rat PTH EIA Sample Diluent – 10 mL Bottle

RUO



Quidel Corporation

2005 State Street, Suite 100

Athens, OH 45701 USA

quidel.com

PI30253100EN00 (11/19)

GLOSSARY

REF

Catalogue number

LOT

Batch code



Use by



Manufacturer



Temperature limitation



Consult e-labeling
instructions for use

RUO

For Research use only

RCNS

Reconstitute with
